**Avaye Danesh Afarin (ADACO)**  
Consulting, Designing & Performing Laboratory Systems

**Avaye Danesh Afarin (ADACO)**  
Consulting, Designing & Performing Laboratory Systems

## ADACO FUME HOOD

### Introduction

Creating and maintaining a safe and healthy environment in the laboratories, which are usually filled with harmful materials, is the vital task expected by each operator working there.

Laboratory fume hood is a type of ventilation system which serves to vent chemical fumes, vapors, gasses, dust, mist and aerosol out of the laboratory and isolate the reaction inside the work zone in order to protect the laboratory from fire and exposures. Also, it enhances safety with creating an air barrier between operator and the harmful materials he/she may handle.

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## Constructions

- Dual wall construction of the fume hood superstructure allows for easy installation of service fixtures.
- Exterior construction built of phosphated steel is resistance to staining and provide a uniformly painted surfaces.
- In spite of improving the cabinet appearance and providing a full coverage, electro static powder coating make the surfaces resistance to abrasion.
- Not collecting contaminants, the cabinet work space has no welded joints. Also there is no screw on the exterior body.
- The superior design of the fume hood with a full accessibility to all the work surfaces provides easy and efficient cleaning.
- Arm rest improves comfort and positions the arms in the level which cannot block the airflow perforations.
- In order to improve the accessibility and visibility of the work chamber, we designed the 10° angled front window.

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## Safety Glass

A frameless sash with Securit glass enhances the safety of the operator and visibility of working area. Securit glass is 5 times more resistant to breakage than normal glass. In case of breakage, there is very low risk for any injury.

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## Microprocessor Control and Alarm System

The microprocessor control is designed to monitor all cabinet functions, display airflow and temperature continuously. The warming-up time, pre-set to ensure the sensors, the fan and the control system are functioning appropriately, can be customized through the control panel. Also, accessibility to the cabinet and the fan settings can be restricted by PIN codes. In case of cabinet dysfunctioning, the alarm system will be immediately aware the operator.

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## Exhaust Blower System

The exhaust blower system is consisted of a direct drive, asynchronous, single or three phase, IP55 motor, a fan motor which is supplied with PPH hubcap and a single block strong high density UV treated and recyclable poly propylene (PPH) housings. The fan performance is conducted in accordance with AMCA 210-85 and IDO 5801.

## OPTIONAL ACCESSORIES/FEATURES

### Worktop

#### Black Natural Granit (standard feature)

Natural granite worktops (thickness of 20-30mm) are chemical resistance, heat proof and easy to clean.

#### Ceramic worktop

Worktop built with ceramic (thickness of 32mm) is anti corrosion, highly resistance to reaction with acids. Moreover, they are solid homogenous materials with no possibility for liquid penetration.

#### Black Epoxy Resin worktop

Epoxy resin is durable, unaffected by most chemicals, flame, heat and moisture, mechanically strong, easy to decontaminate and above all will not delaminate, swell or crack.

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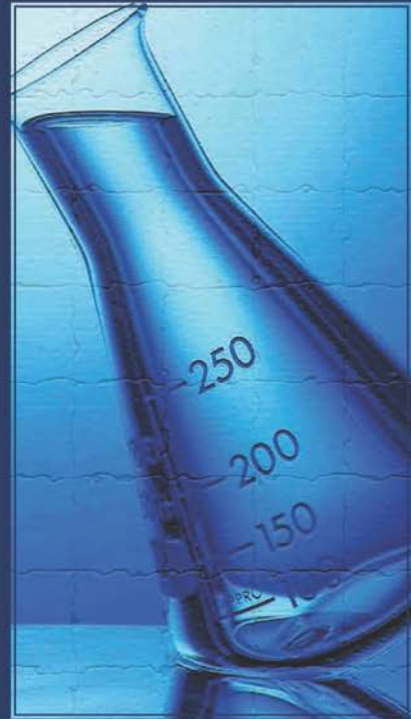
### Cup Sinks

#### a. Polyester Resin Cup Sink(standard feature)

Installed in size of 300×150 mm, poly cup sinks are chemical resistant, easy to clean, unaffected by acids, alkalies and boiling solutions.

#### b Ceramic Cup Sink

Available in size of 300×150 mm, ceramic cup sinks are non-porous, heat resistance and won't rust or discolor.

### Laminated Safety Sash Glass

Manufacturing with an interlayer polyvinyl butyral (PVB) between two layers of glass, the laminated glass is held in place when breaking and does not break to sharp pieces. Thus, operator safety is guaranteed.

### Fittings

The standard cabinet comes with one remote control water and one remote control gas tap. The operator can enjoy working while the nozzles are located inside the work chamber. Compressed air and vacuum taps can be added in case of customer demand.

### 8. Fixtures

type		type		type		type		type		type	
A		B		C		D		E		F	
G		H		I		J		K			

### Distillation Grid

Maximizing usable work area, glassware and other small apparatus can be mounted to the rear wall of work zone through distillation grid built with stainless rods.

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### Base Cabinet

The base cabinet is available in phosphate steel with iron profiles coated by electrostatic powder. The cabinet is equipped with adjustable legs. For easy plumbing installation, an opening is designed on the back wall. In addition, the operator can enjoy a big leg room as the base cabinet depth is 100 mm less. Customer can select a cabinet with support stand and removable base cabinet.

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## FUME HOOD SELECTION

To select a fume hood, following recommendations should be considered:

### 1. The type of operation in the hood

- If perchloric acid or radioactive materials are handled, special fume hood should be ordered.
- Large apparatus or complex set-ups may require big work zone. So, walk-in fume hoods or low bench fume hoods are the recommended options.
- In case of working with large volume of flammable substances, the hood should be equipped with explosion-proof lights, a fire-supervision system and ATEX exhaust fan.

### 2. Air consumption of hood

Depend on the laboratory conditions, you are recommended to choose special type of hood:

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Type of hood	Conditions	
	The room air is adequate	The outside temperature is extreme
Bypass Hood	Yes	No
Variable Air Volume Hood	Yes/No	Yes
Add Air Hood	No	No

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## FUME HOOD OPERATION

1. When you stand in front of a laboratory fume hood, a turbulent area is formed in front of you. Thus, it's seriously recommended to keep materials at least 6 inch inside the hood.
2. The protection is greater as much as you place the materials farther behind the sash.
3. The large equipments should be placed on the legs not to cause dead spots or reverse airflow.
4. Do not cause airflow disturbance by making rapid movements inside and outside the hood or walk quickly past the hood.
5. The sash should be closed if you don't need to access inside the hood continuously.
6. The hood is not a storage cabinet. So, keep the unusable materials outside the work zone.
7. Personal protective equipments like splash goggles or gloves should be wearing when operating.

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## FUME HOOD LOCATION

As the performance of a fume hood is greatly affected by the surrounding conditions, the hood should be located away from open doors, windows, the line of traffic and high velocity air diffusers.

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