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

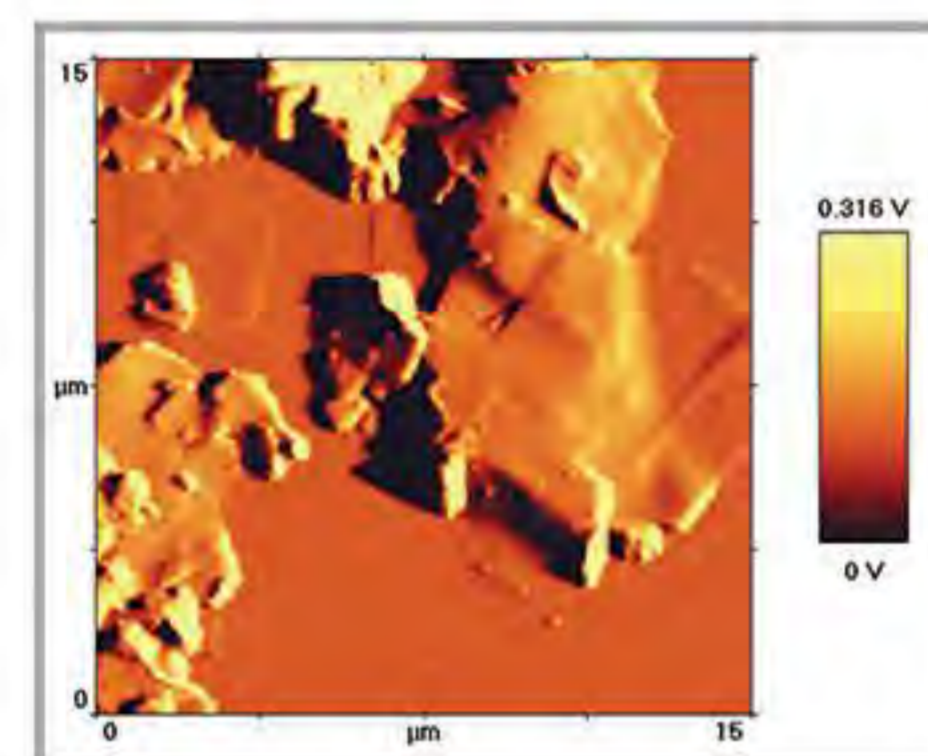
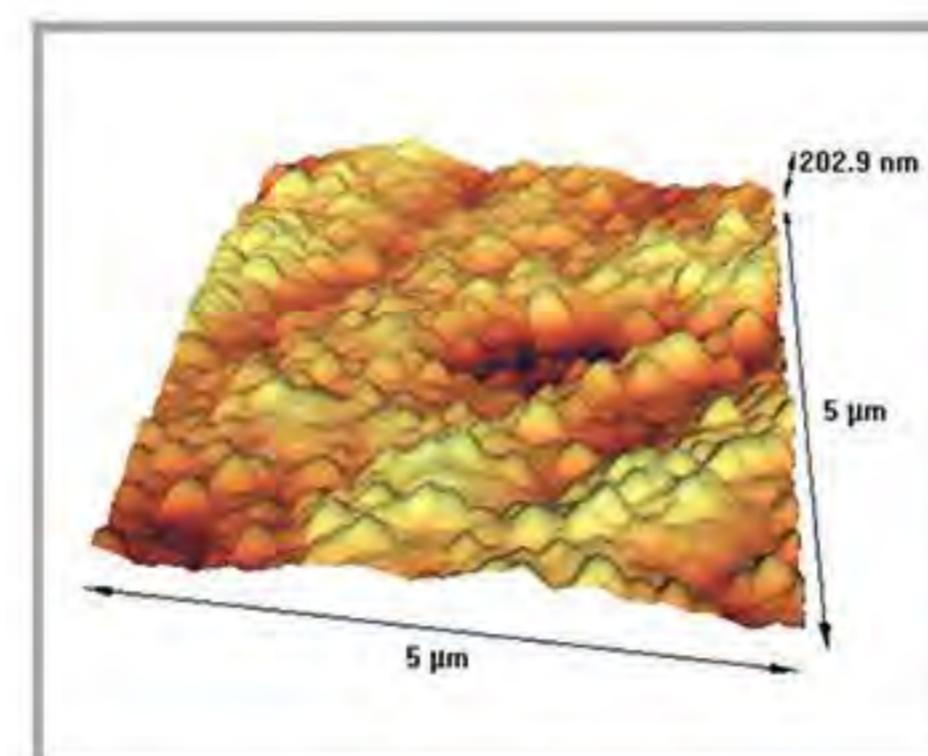
Gold nano-sheets
Dr. H. Dehdashti
University of Shiraz

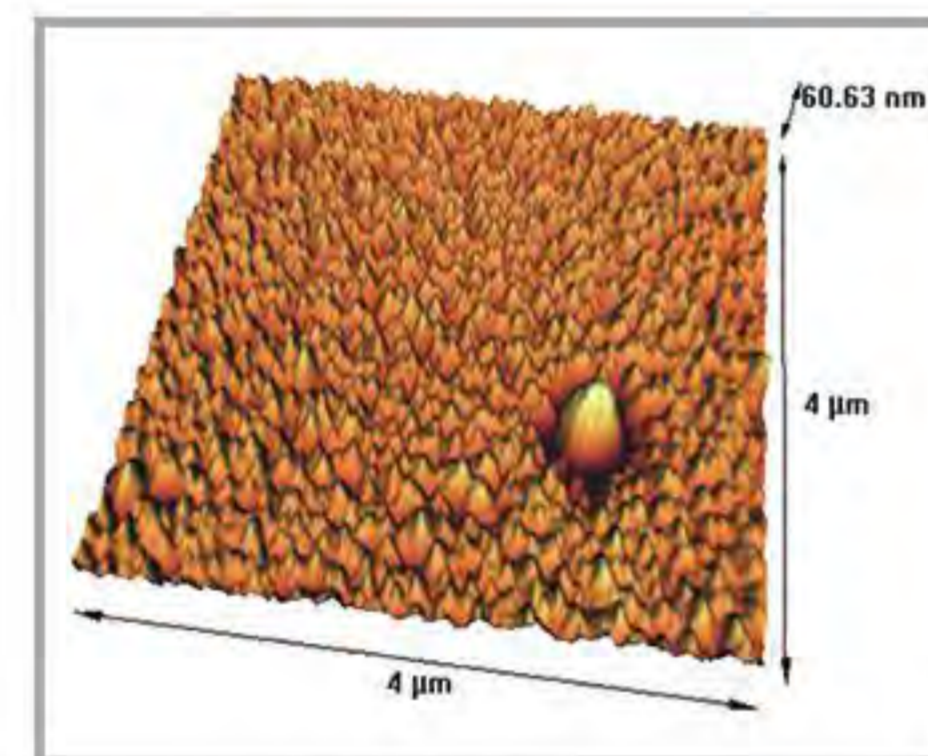
Iron nitrates coating on Al by Active Screen Plasma Nitriding (ASPN) Method
Dr. M. Soltaniyeh, IUST

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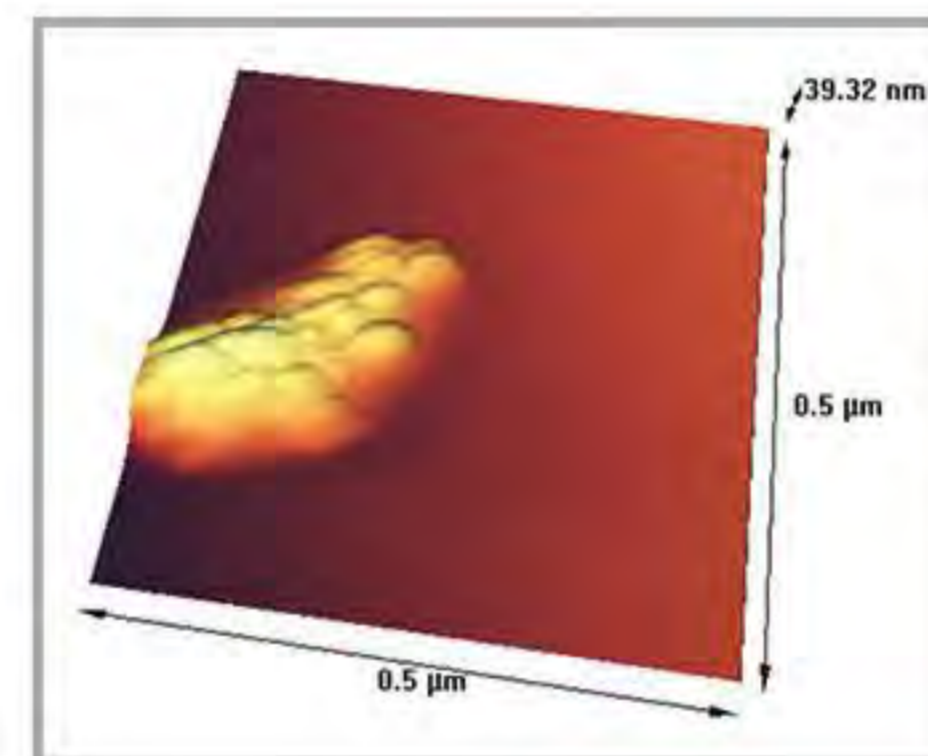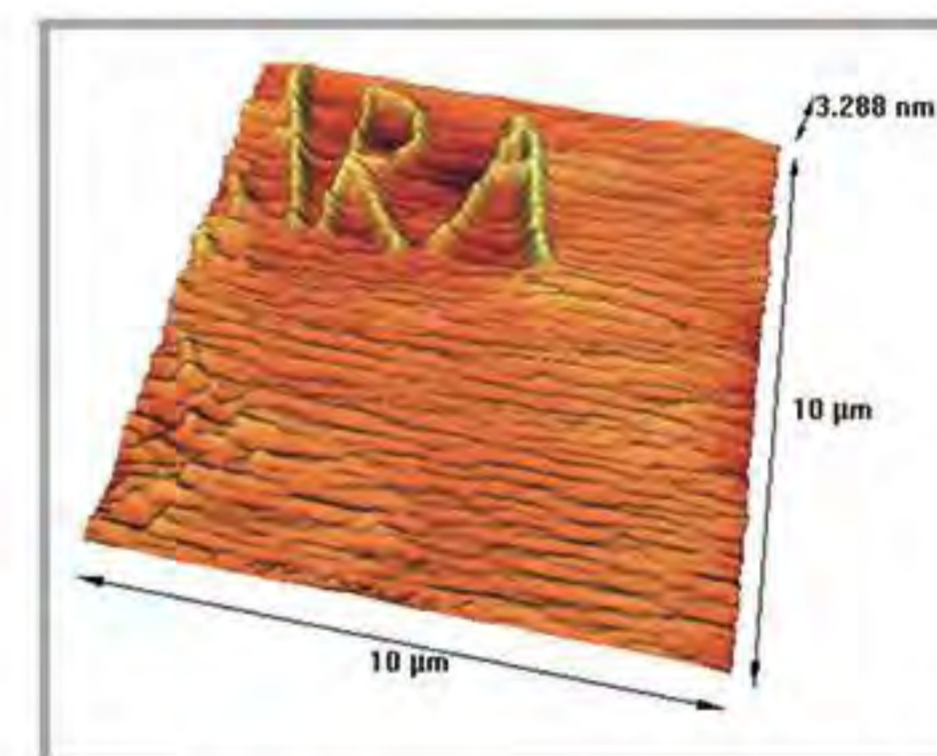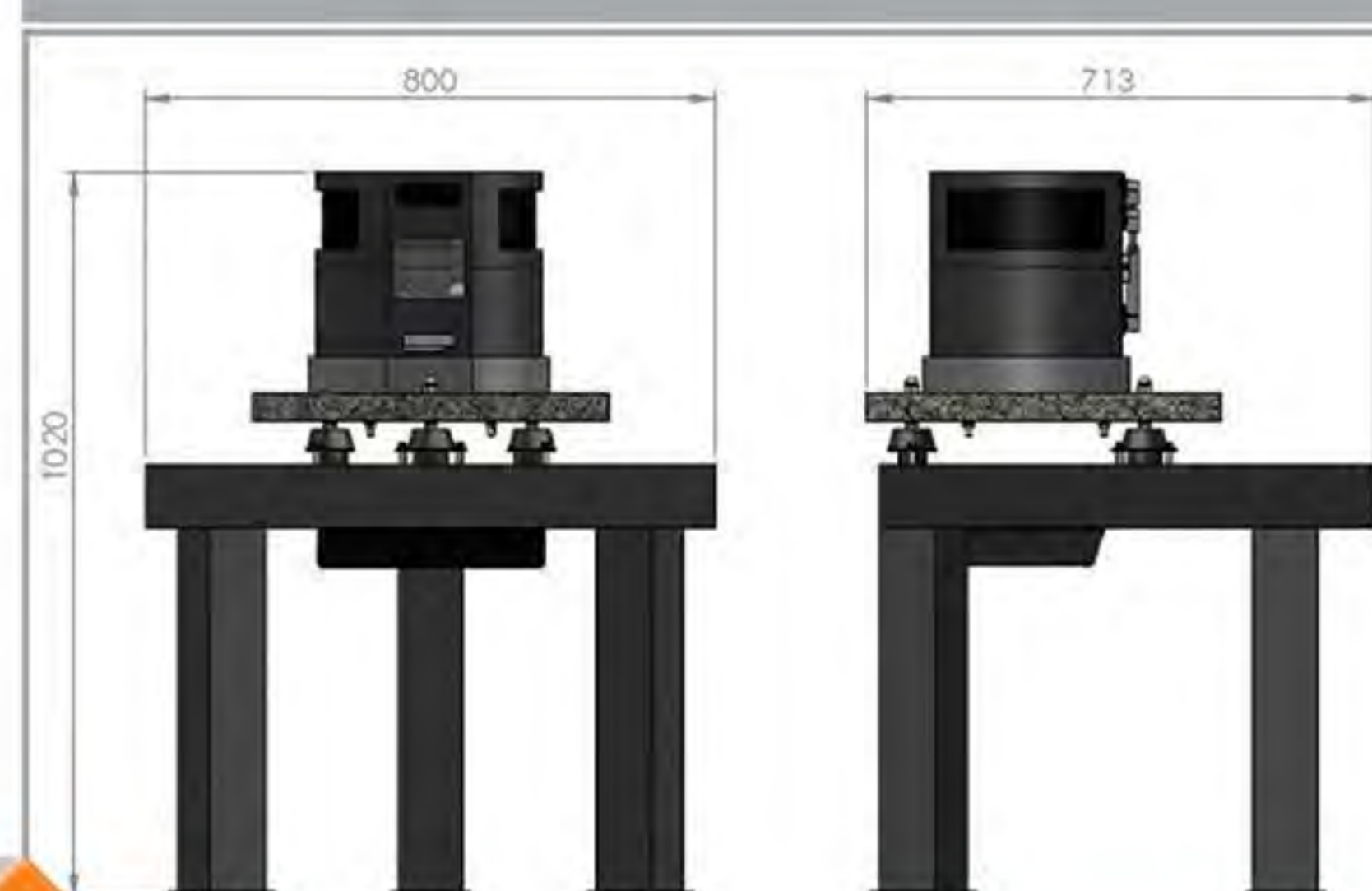
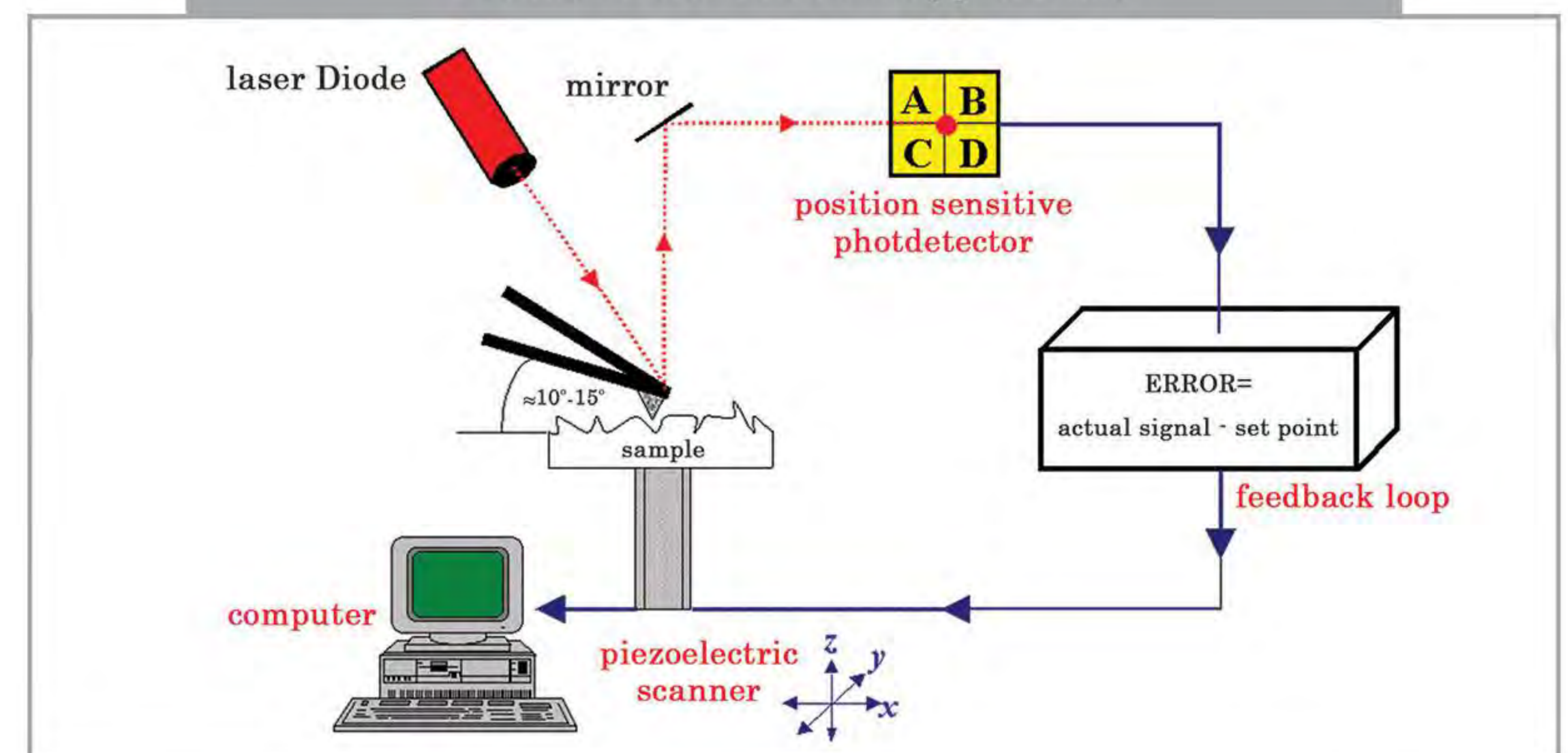
- Contact Mode
- Dynamic Mode
- Tapping Mode
- Frequency Modulation

Nano Properties Determination

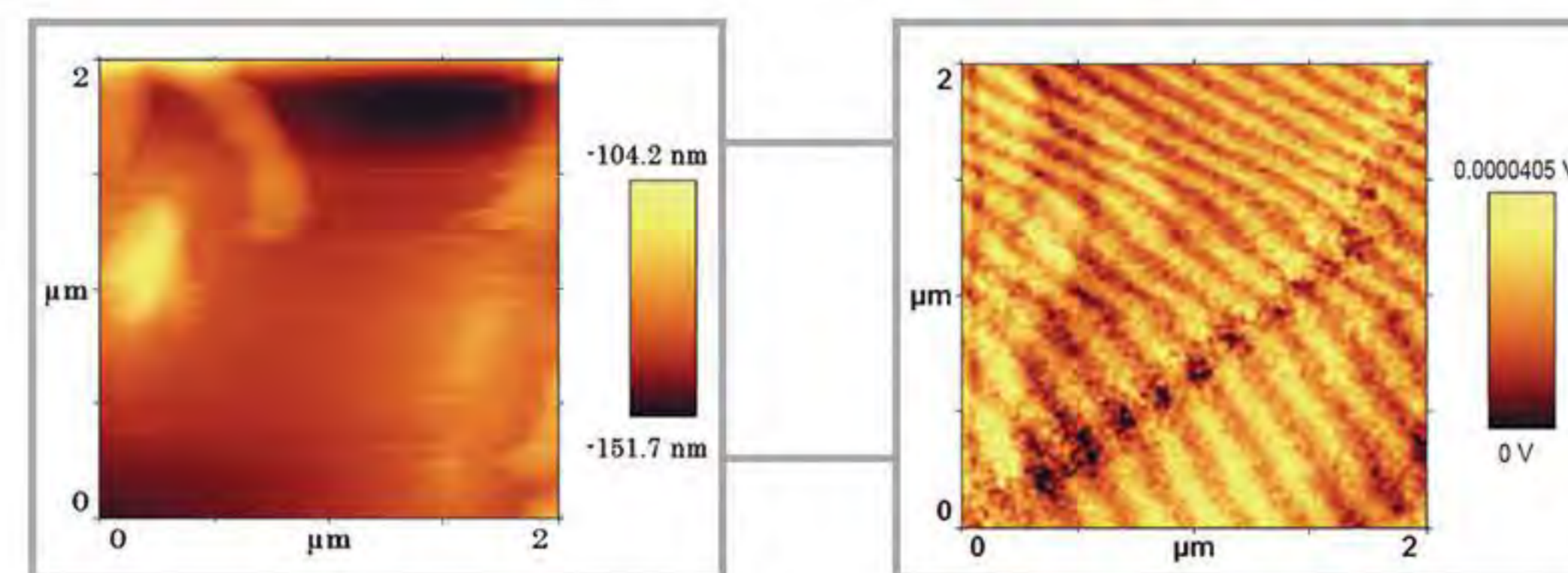
- Magnetic Force Microscopy (MFM)
- Electric Force Microscopy (EFM)
- Conductive AFM (CAFM)
- Kelvin Probe Force Microscopy (KPFM)
- Piezoresponse Force Microscopy (PFM)
- Lateral Force Microscopy (LFM)
- Force Modulation Microscopy (FMM)
- Force Spectroscopy

Nano Lithography and Manipulation


- Chemical Nano-Lithography
- Mechanical Nano-Lithography

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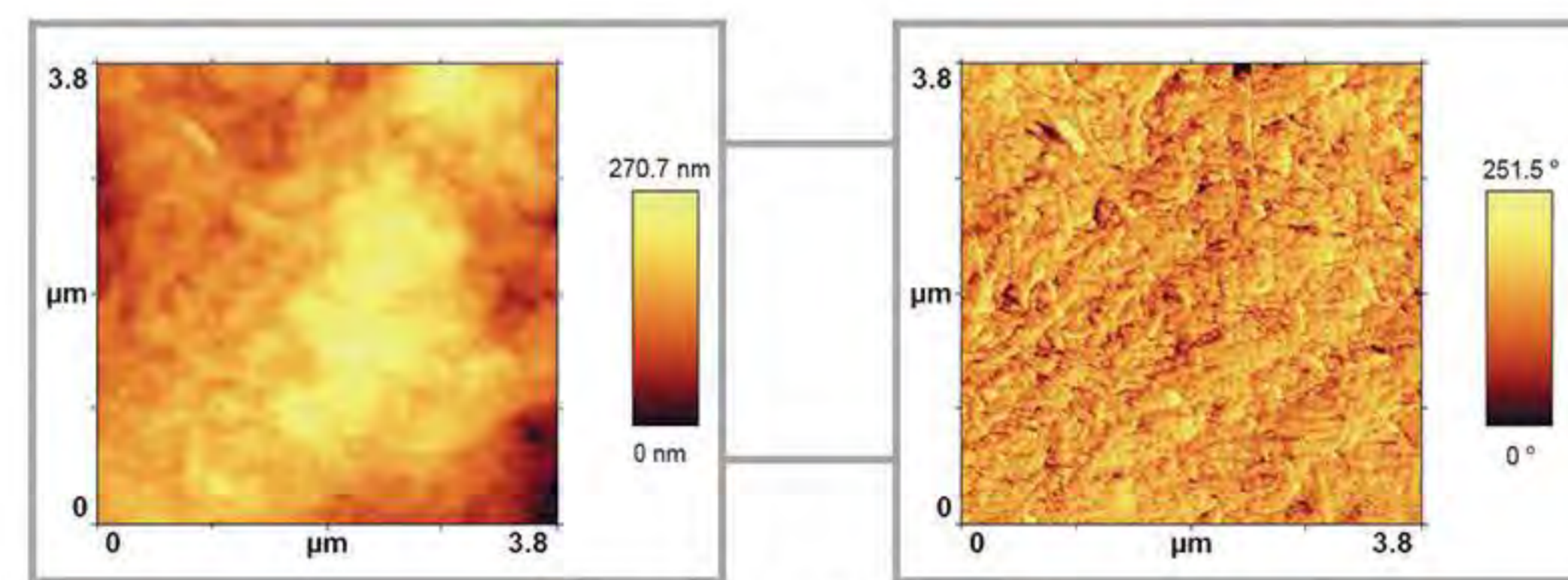
Topo (left) and MFM (right) images of FeSiB Magnetic Ribbon
Dr. M. Soltanieh, IUST

Phase images of human tooth
Dr. Y. Seyed Jalili
Islamic Azad University Science and Research Branch

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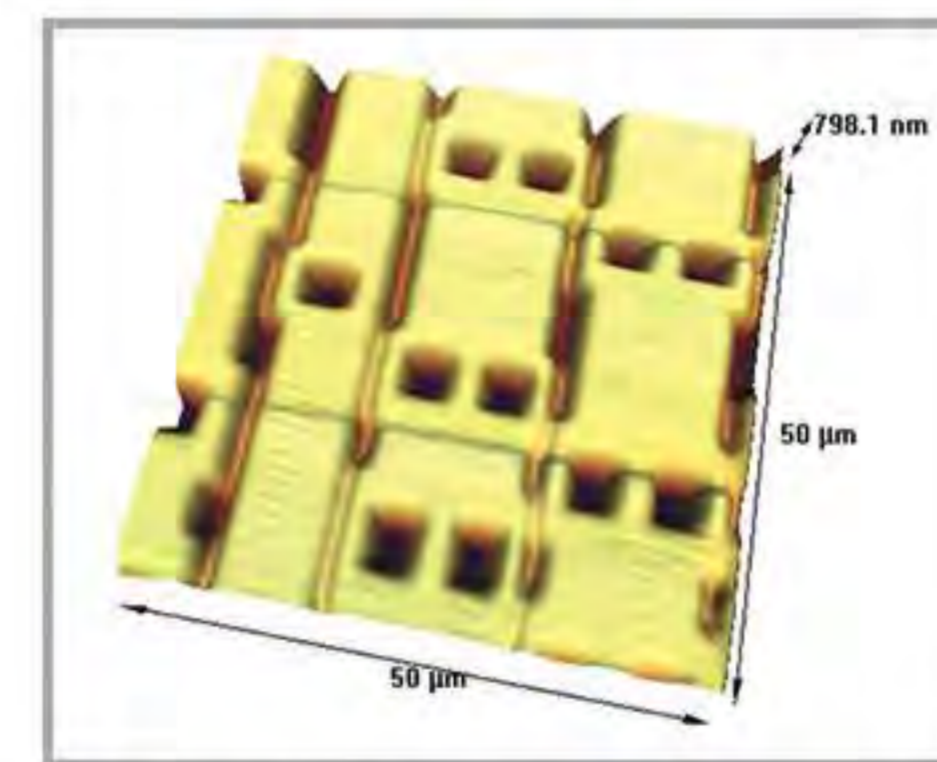
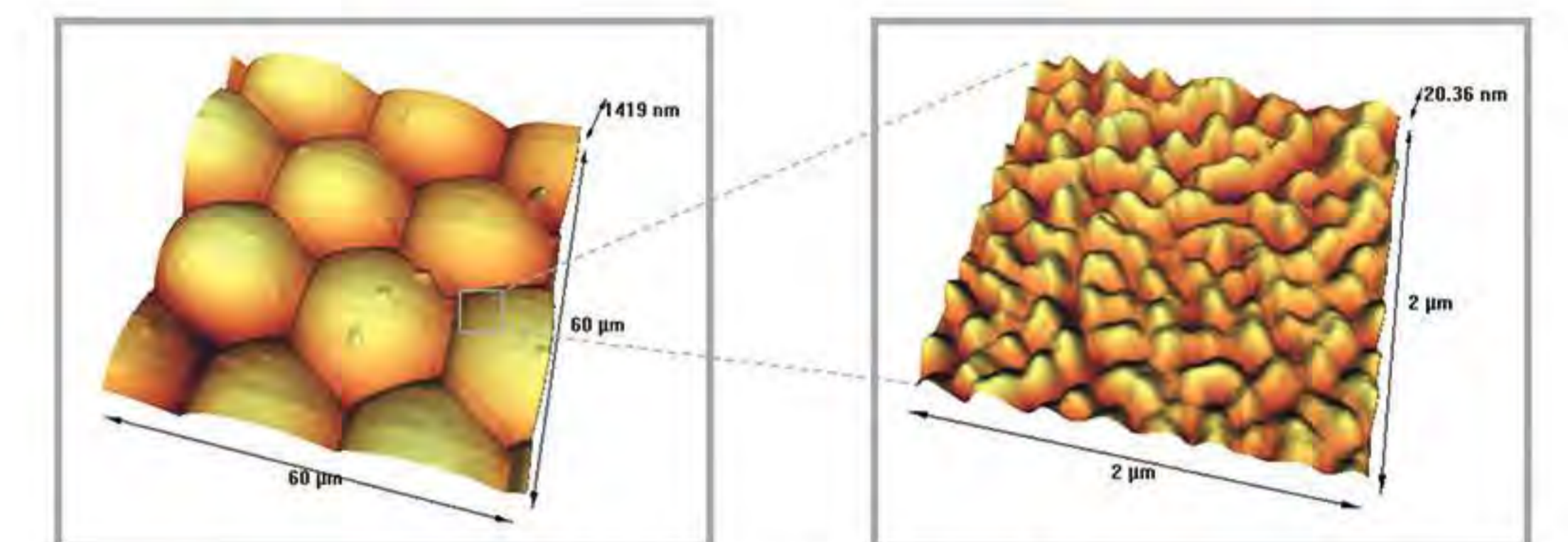
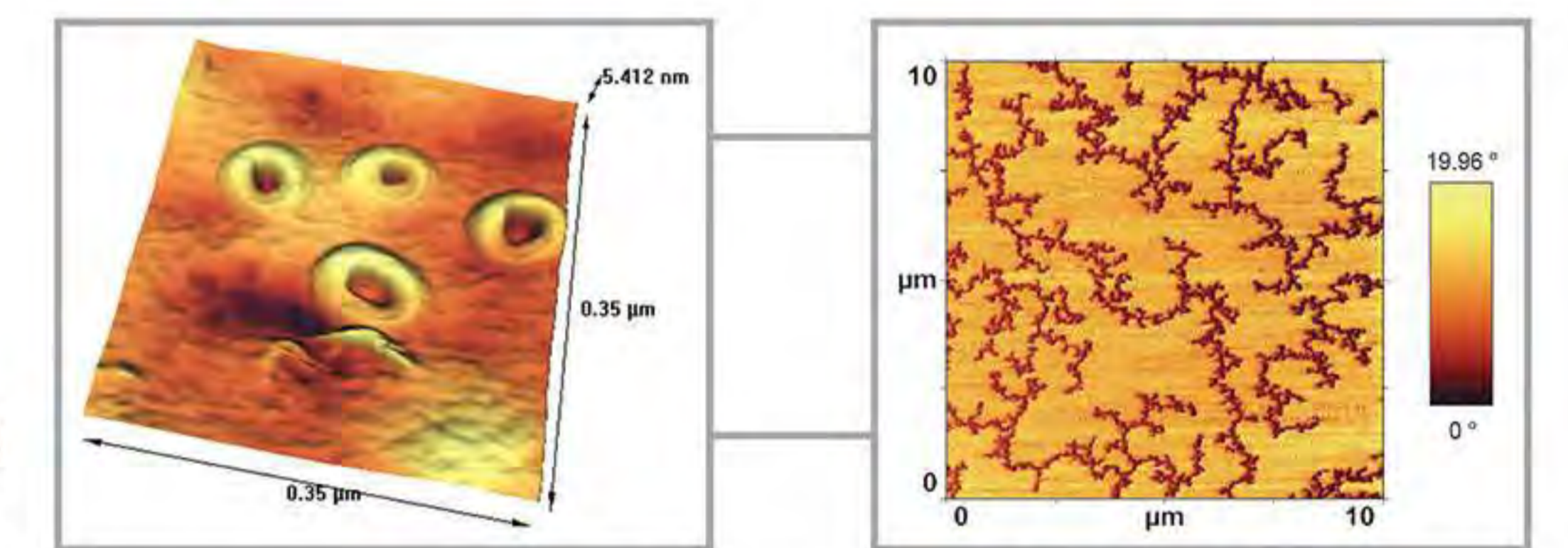
Bee's eye, Mr. H. Akbari, Ara Research Centre

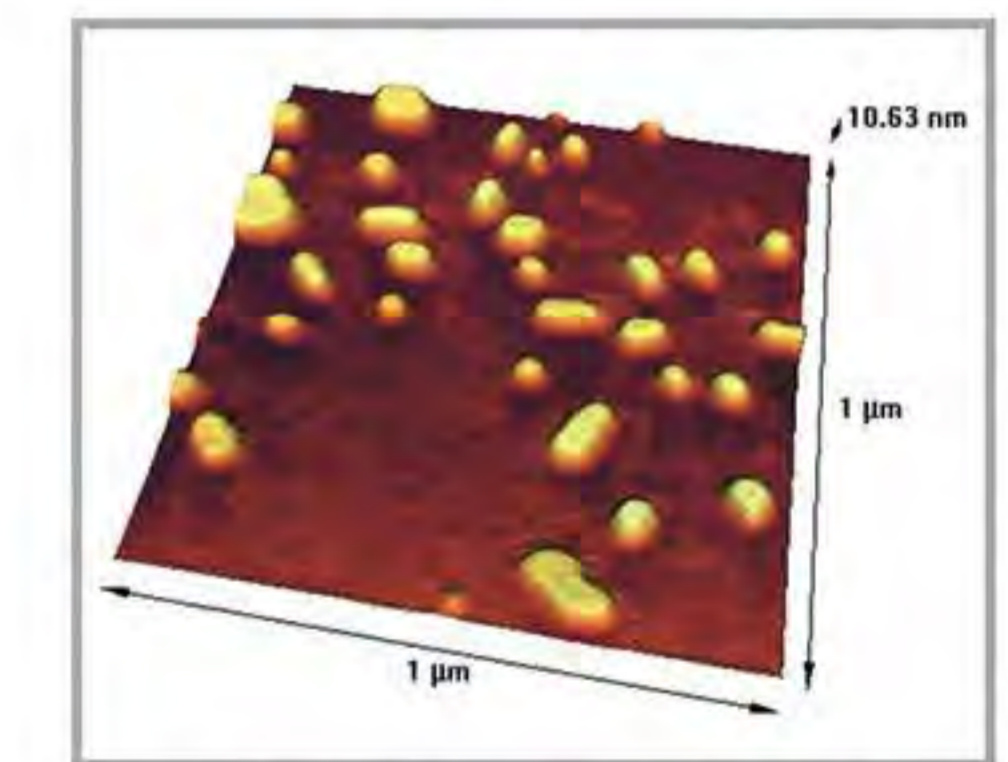Tau Protein + ps-9 (left) and Tau Protein + pn-8 (right)
Dr. M.R. Lornejad
Anahita Life Science

Influenza virus, Dr. E. Saber Far
Ara Research Centre

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	<table border="1"> <tr><td>Head</td><td>H-A11</td></tr> <tr><td>Scanner</td><td>S-C2</td></tr> <tr><td>Controller</td><td>C-2NFP</td></tr> <tr><td>Options</td><td>HighSpeed Module, Motorized Head H-A11 S-C3, S-C4, S-C5</td></tr> <tr><td>Vacuum Pump</td><td>Compatible with well-known pump e.g. Leybold TriVac</td></tr> <tr><td>Vacuum Gauge</td><td>Compatible with well-known gauge e.g. Leybold Pirani gauge</td></tr> </table>	Head	H-A11	Scanner	S-C2	Controller	C-2NFP	Options	HighSpeed Module, Motorized Head H-A11 S-C3, S-C4, S-C5	Vacuum Pump	Compatible with well-known pump e.g. Leybold TriVac	Vacuum Gauge	Compatible with well-known gauge e.g. Leybold Pirani gauge
Head	H-A11												
Scanner	S-C2												
Controller	C-2NFP												
Options	HighSpeed Module, Motorized Head H-A11 S-C3, S-C4, S-C5												
Vacuum Pump	Compatible with well-known pump e.g. Leybold TriVac												
Vacuum Gauge	Compatible with well-known gauge e.g. Leybold Pirani gauge												

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
	<table border="1"> <tr><td>Head</td><td>H-M21</td></tr> <tr><td>Scanner</td><td>S-NC2</td></tr> <tr><td>Controller</td><td>C-2BFP</td></tr> <tr><td>Options</td><td>HighSpeed Module</td></tr> <tr><td>Optical Microscope</td><td>Included Inverted Microscope</td></tr> </table>	Head	H-M21	Scanner	S-NC2	Controller	C-2BFP	Options	HighSpeed Module	Optical Microscope	Included Inverted Microscope
Head	H-M21										
Scanner	S-NC2										
Controller	C-2BFP										
Options	HighSpeed Module										
Optical Microscope	Included Inverted Microscope										

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
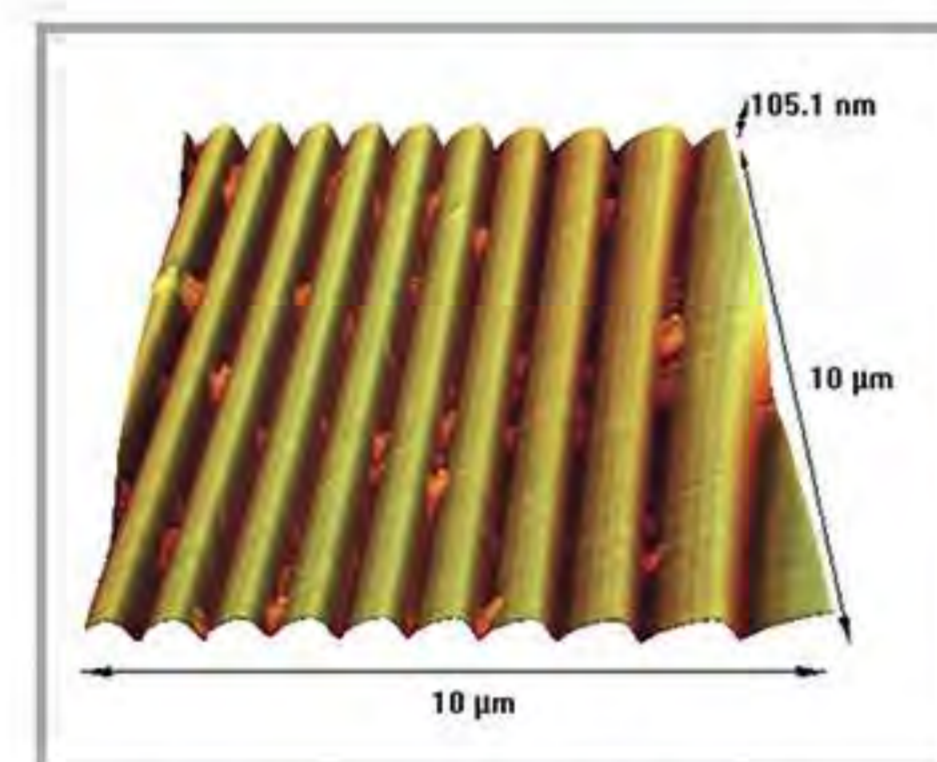
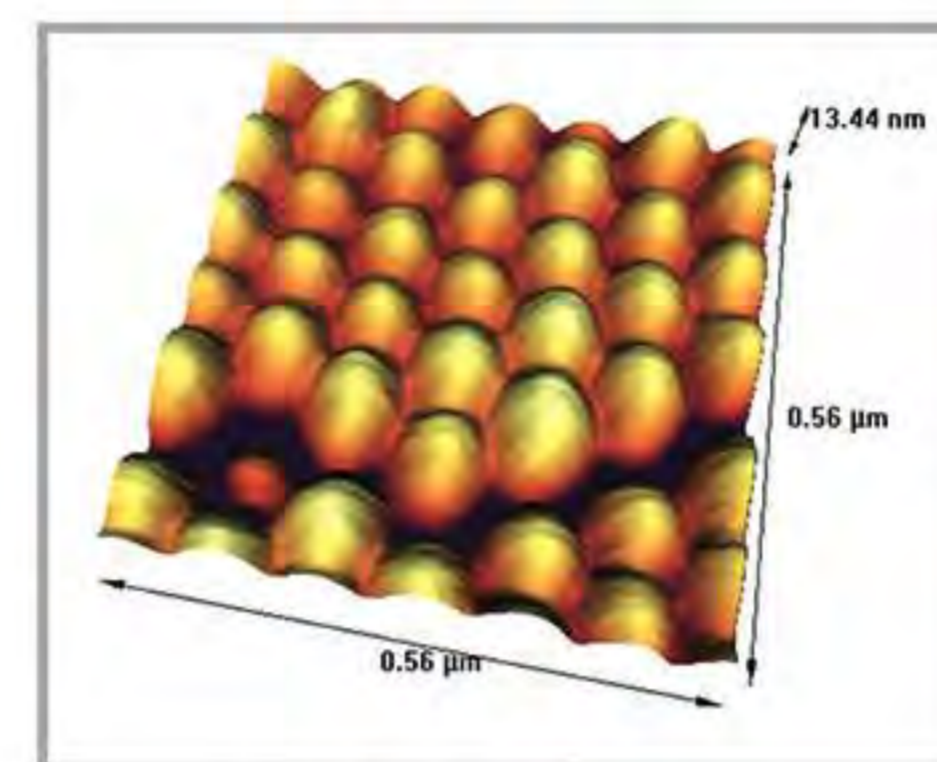
Ag coating on DVD, Dr. Hamidi
Shahid Beheshti University

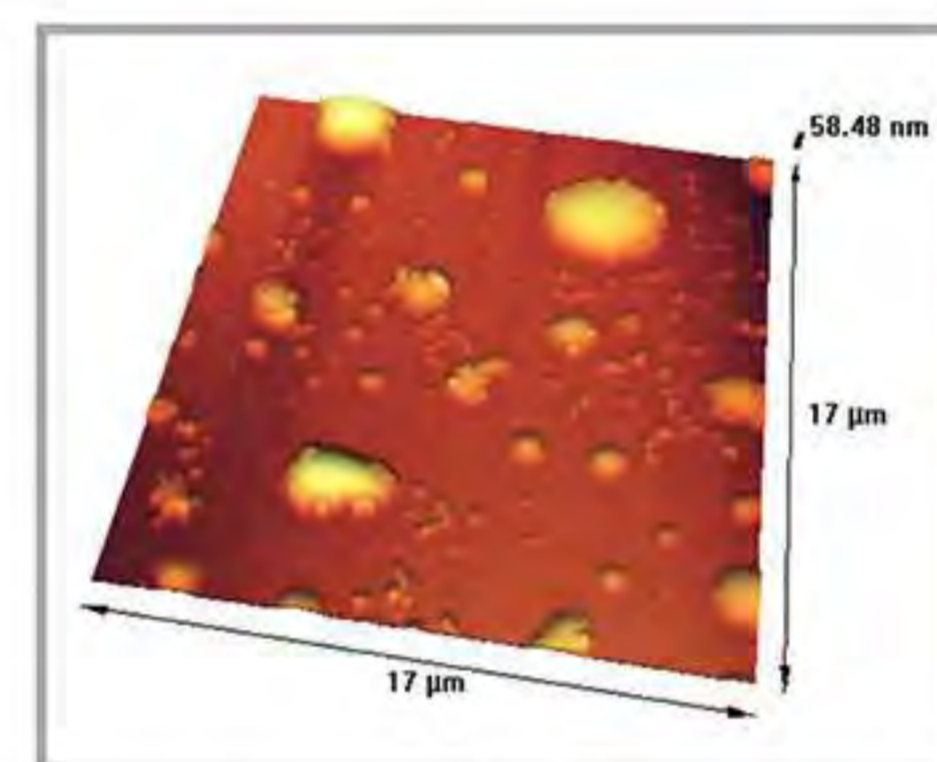
Latex, Ms. Z. Sarkhosh
Ara Research Centre

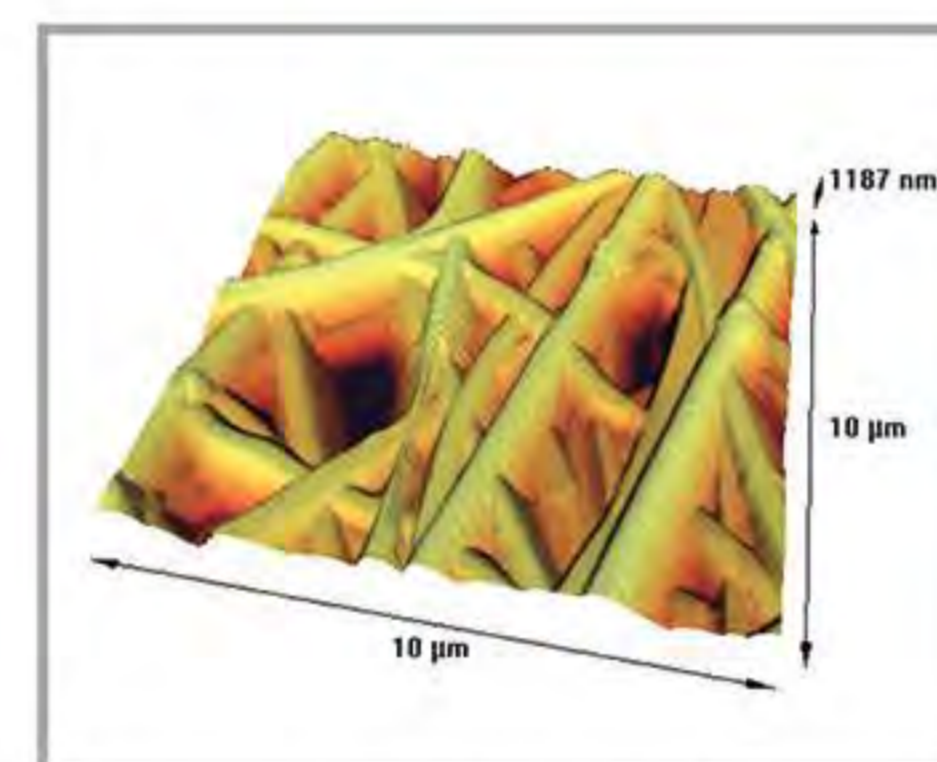Liposome, Ms. F. Khakrah
Ferdowsi University of Mashhad

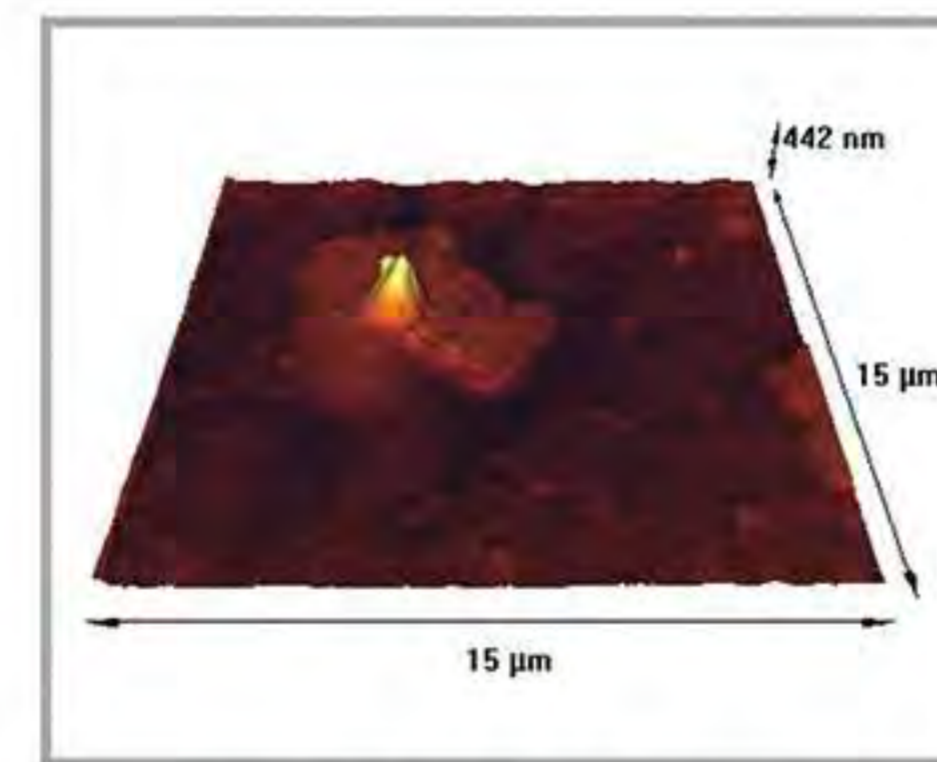
Electrospun fibers, Dr. M. Ghomi
Islamic Azad University
Pharmaceutical Sciences Branch

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
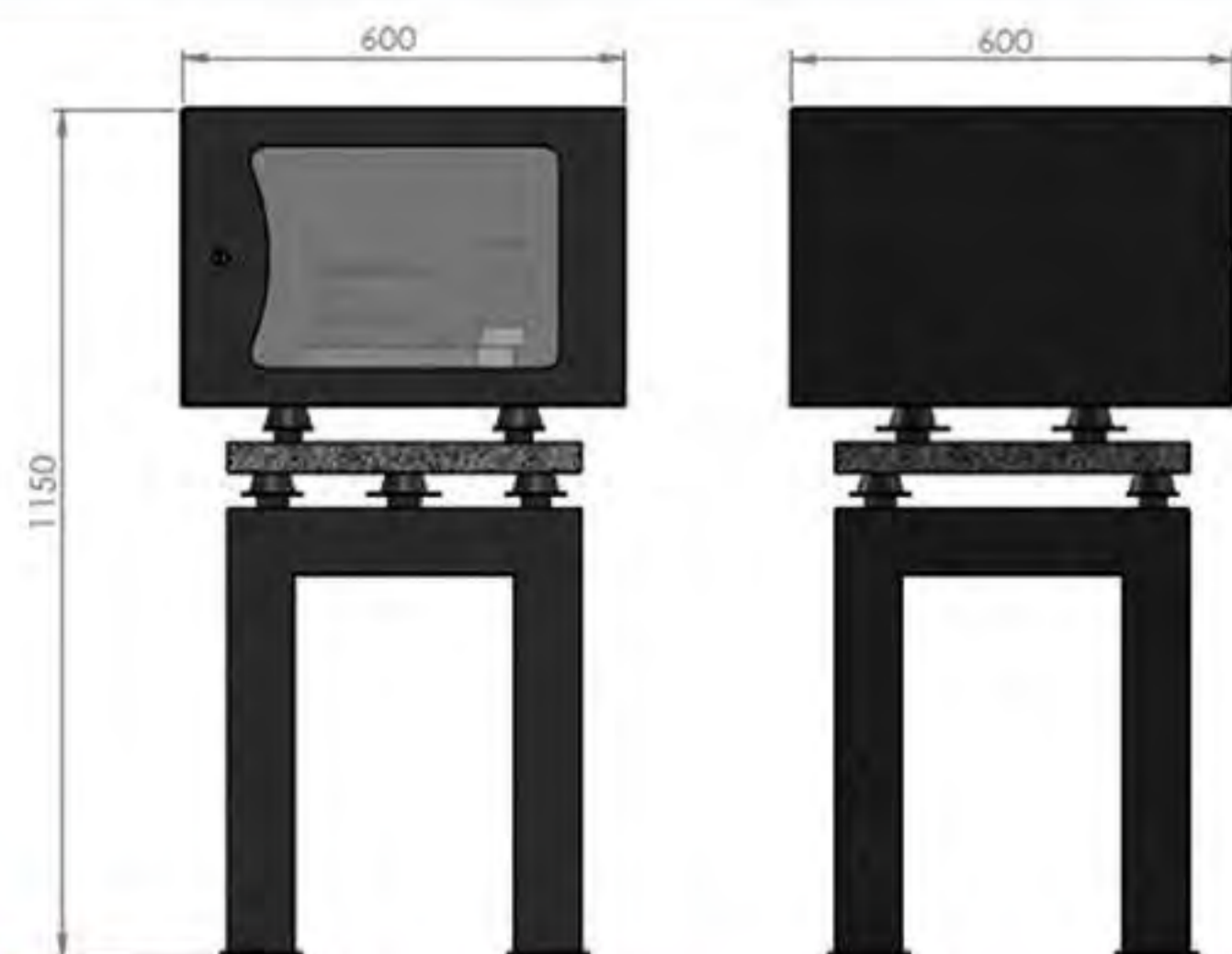
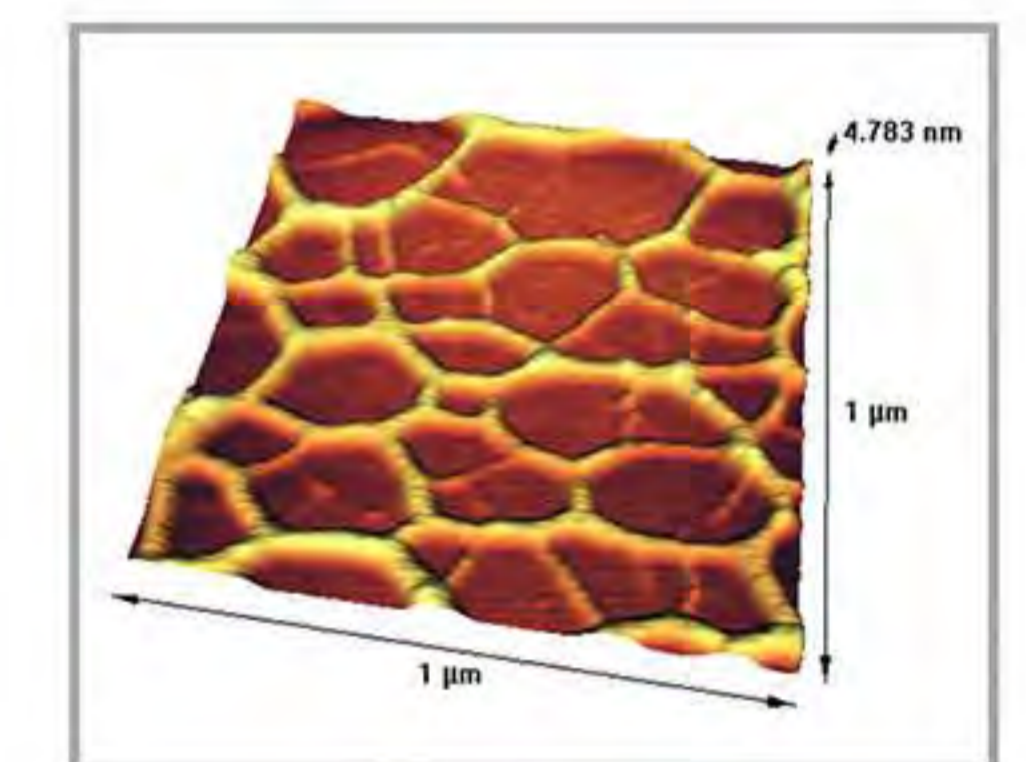
Albumin protein
Ms. F. Khakrah
Ferdowsi University of Mashhad

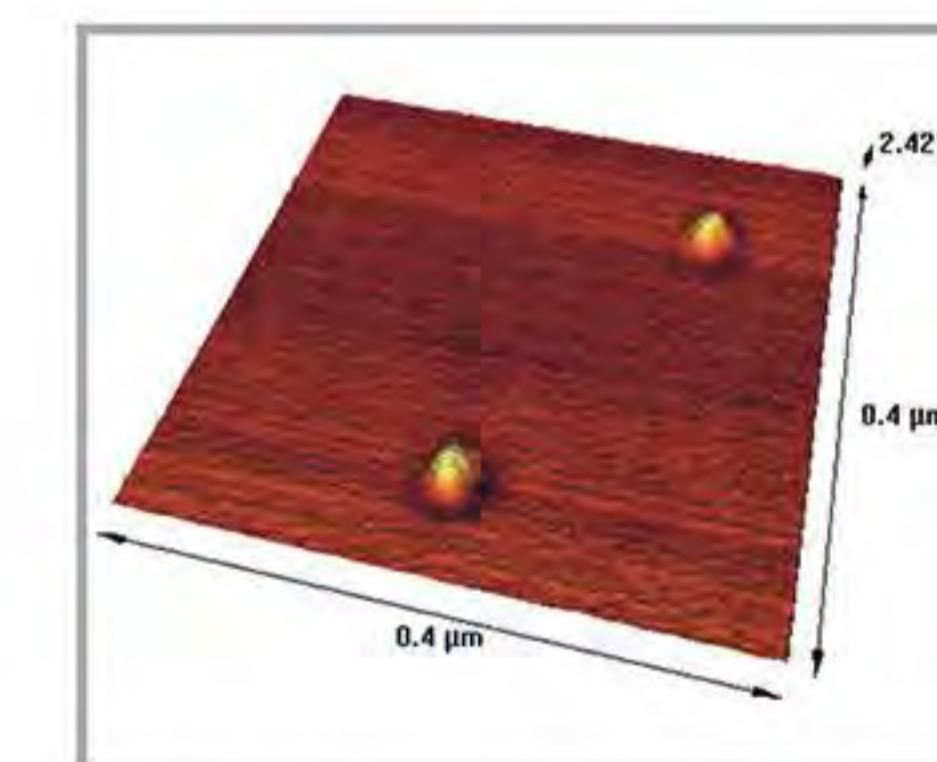
DNA network
Dr. M.R. Lornejad
Anahita Life Science

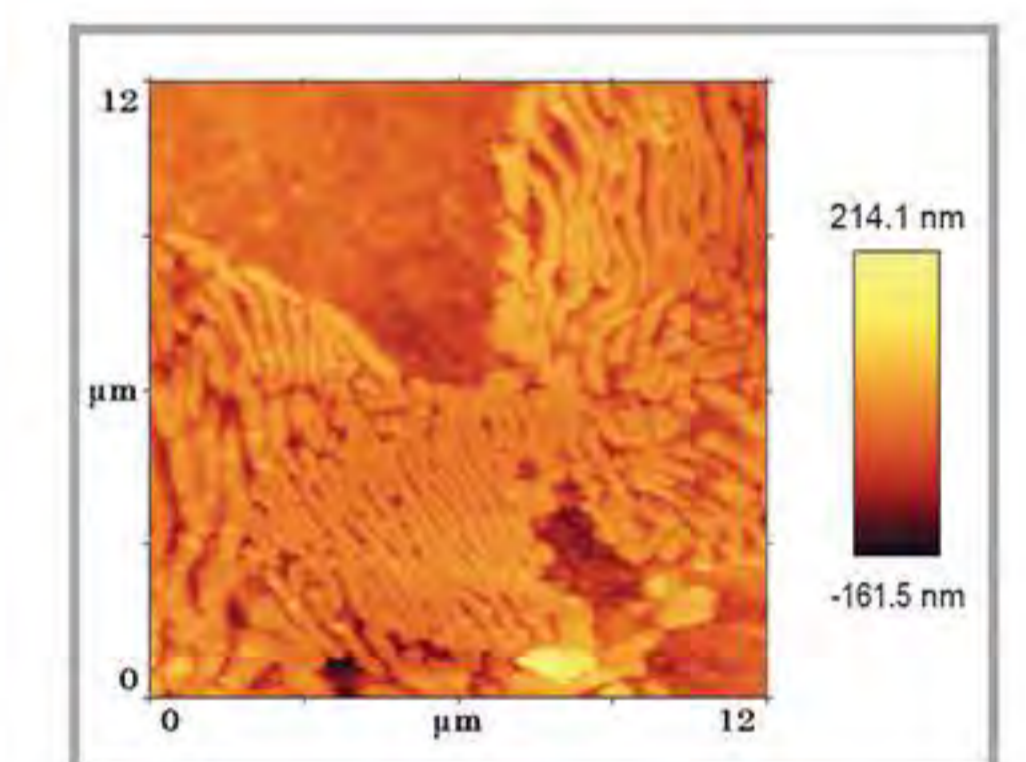ZnO + Cr nano-particles phase image
Ms. S. Kerdegari
Ara Research Centre

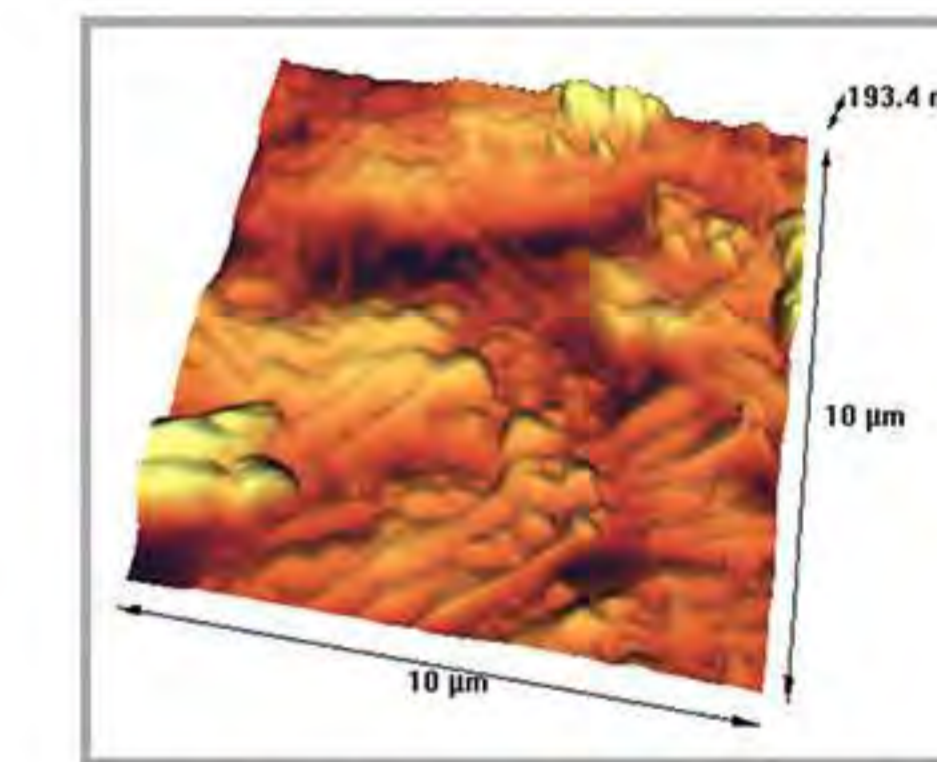
Metallic alloy micro structure
(Pearlitic SPK)
Mr. H. Akbari, Ara Research Centre

Rabbit cartilage, Ms. F. Khakrah
Ferdowsi University of Mashhad

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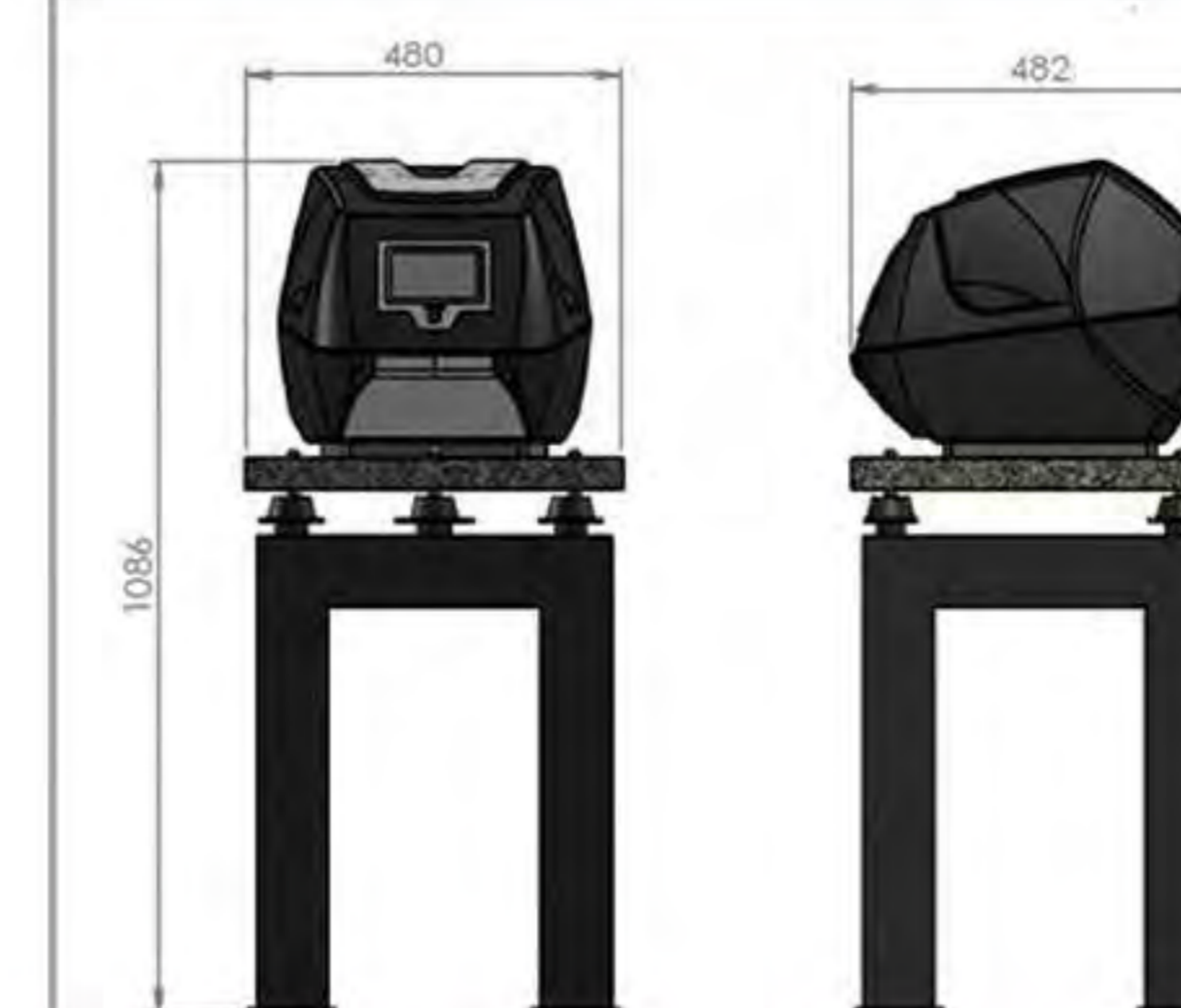Head	H-M11
Scanner	S-C2
Controller	C-2HFP
Options	Motorized Head H-A11 S-C3, S-C4, S-C5

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Head:

- Laser diode maximum 3 mW, 670 nm with lens system
- Integrated 4-quadrant photo-detector with amplifier electronics
- Laser beam adjustment on the photo detector in two directions
- Built-in dither piezo for acoustical excitation in dynamic mode
- Color-camera with microscope optic with a direct view onto the cantilever: 640x480 pixels (VGA)

Model	H-M11	H-A11	H-M21
Optical Adjustment	Manual	Auto	Manual
Approach method	Auto/Manual	Auto/Manual	Auto/Manual
Top view camera	20X up to 200X zoom	20X up to 200X zoom	20X up to 200X zoom

Scanner:

Scanners of various x-y range and z displacement are produced by Ara Research. Scanners features are listed in the following table

Model	S-C1	S-C2/S-NC2	S-NC3	S-C4	S-C5
Scan range XY	30 μm	50 μm	80 μm	50 μm	80 μm
Scan Range Z	3.5 μm	3.5 μm	3.5 μm	7 μm	7 μm
Resolution XY	2 nm	1 nm	1 nm	1 nm	1 nm
Resolution Z	0.3 nm	0.1 nm	0.1 nm	0.1 nm	0.1 nm
XY sample movement	6.5 mm	6.5 mm	6.5 mm	6.5 mm	6.5 mm
Max Sample Thickness	7 mm	7 mm	7 mm	7 mm	7 mm

Controller:

2nd generation of ARA-AFM controller improved in:

- 1- Size reduction
- 2- Developing control capability
- 3- Noise reduction and stability improvement

Controller model No. is detailed as follows:

Device Model	Supported Modes
E ducational	S tandard
M ulti model	A dvanced
B io	F ull
N anoVac	F ull P lus
H ighSpeed	

Software:

Number of scanning channels	up to 8
Number of points per image	up to 1024x1024
Number of points per spectrum	up to 1024
Angular scanning	Available
Resume scan	Available
Oxidation lithography	Available
Multi-scan	Available
Operating System	Microsoft Windows XP, Windows 7, Windows 8, Windows 10
Program interface (SDK)	Nano Ara & Imager

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