

## شرایط محیطی مورد نیاز برای نصب دستگاه Ethidium Bromide Removal

### Instalation Instructions

1. Place the Extractor device on a sturdy and level surface near the sink.
2. Put the drain tube in a storage container or in the sink
3. Upon completion of filtration, note the amount of as a record of the total amount of buffer decontaminated with each cartridge.
4. The filtrate can be safely discarded by pouring down the sink, since up to 99% of the EtBr has been adsorbed by the Extractor device.
5. After 1000 liters of staining buffer have been decontaminated, dispose of the Extractor cartridge in accordance with regulations set by the institution or locality for solid hazardous waste disposal.

### Maintenance

The ET Adsorb may be cleaned with a moist cloth containing a mild soap solution. The blocks are aluminum and may be cleaned with any of the commercial aluminum cleaners on the market.